

1. Search for the gene(s) of interest on The RNAi Consortium (TRC) portal (<http://www.broadinstitute.org/rnai/public/>) and select the ~five highest scoring non-overlapping targeting sequences.
 - *In general, sequences targeting the CDS are preferred because they can be rapidly screened by transient co-transfection with an expression plasmid for the gene to be knocked down*
 - *Sequences that are 100% human specific are preferred so that future addback experiments can be performed with the mouse or rat homolog*
 - *Custom targeting sequences for special applications can be designed according to the TRC design rules: <http://www.broadinstitute.org/science/projects/rnai-consortium/trc-shrna-design-process>*
2. Order the primers (minimum synthesis, standard desalting) listed under “Oligo design for arrayed cloning” after replacing the XhoI site in the loop (CTCGAG) with a PstI site (CTGCAG). Dissolve primers to a concentration of 50 μ M in PCR-grade H₂O
 - *Unlike XhoI, PstI restriction sites are absent in pLKO.1, allowing ligated hairpins to be screened rapidly by restriction digest*
3. Anneal the shRNA primers by setting up the following reaction in a PCR tube (one annealing reaction per shRNA duplex):
 - 2.5 μ l top-strand primer
 - 2.5 μ l bottom-strand primer
 - 5 μ l 10 \times annealing buffer
 - 40 μ l ddH₂O
 -
 - 50 μ l total volume
4. Heat the annealing reaction to 95°C for 5 min in a PCR thermocycler. Unplug the thermocycler and allow the primers to anneal by cooling to room temperature overnight.
 - *The shRNA oligos are designed to contain restriction enzyme-like cleavage ends after annealing; therefore, they do not need to be digested*
5. Phosphorylate the annealed shRNA primers by setting up the following reaction in a PCR tube:
 - 1 μ l annealed primers
 - 1 μ l 10 mM ATP
 - 1 μ l 10 \times PNK reaction buffer (supplied with NEB #M0201)
 - 1 μ l T4 polynucleotide kinase (NEB #M0201)
 - 6 μ l ddH₂O
 -
 - 10 μ l total volume
6. Incubate the phosphorylation reaction at 37°C for 30 min, then heat inactivate at 70°C for 10 min.
7. In parallel with the oligo annealing and phosphorylation, set up three 50 μ l restriction digests, each containing 2 μ g pLKO.1 or tet-pLKO.1, 5 μ l 10 \times EcoRI buffer, 2.5 μ l EcoRI, 2.5 μ l AgeI, and 0.5 μ l of 100 \times BSA.
 - *We have DNA stocks of pLKO.1 puro, pLKO.1 neo, pLKO.1 hygro, and tet-pLKO.1 puro which can be used depending on the combination of perturbations desired*
8. Digest pLKO.1 or tet-pLKO.1 at 37°C for 3 hr and purify the ~7 kb (if pLKO.1) or ~9 kb (if tet-pLKO.1) digested fragment on a 1% agarose gel
 - *The pLKO.1 series releases a 1.9 kb “stuffer” sequence after EcoRI-AgeI digest, which readily separates the double-cut vector from any single-cut vector*
9. Excise the digested vector, purify, and ethanol precipitate the three digests into 5 μ l EB as described in Janes_PCRcloning.pdf.
10. Set up a 10 μ l ligation reaction containing 7.5 μ l phosphorylated-annealed shRNA oligo, 100–250 ng digested pLKO.1, 1 μ l 10 \times T4 ligation buffer, and 0.5 μ l T4 DNA ligase (New England Biolabs #M0202S).
 - *Be sure that the ligation buffer has thawed completely (no visible precipitates) and is kept at room temperature before use*
 - *Note this reaction volume is half that recommended by the manufacturer*

shRNA cloning into pLKO.1 vectors

Janes Lab Protocols

Entered by Karin Holmberg, Kevin Janes

12/14/14

- *Be sure to include a control ligation that includes everything but the shRNA oligo*
11. Mix by pipetting gently, incubate at room temperature for 1 hr, and then ethanol precipitate and transform into electrocompetent bacteria as described in Janes_PCRcloning.pdf.
 - *tet-pLKO.1 cultures require incubation at 30°C to minimize recombination of the plasmid*
 12. Perform a PstI-BamHI diagnostic digest on ~four ampicillin-resistant colonies per shRNA oligo. Positive clones will yield two products at ~6.3 kb and ~700 bp (for pLKO.1 vectors) and ~6.2 kb and ~2.6 kb (for tet-pLKO.1).
 13. Send three digest-positive clones out for forward shRNA sequencing at Genewiz with the “hairpin” option on the website.
 14. Screen sequence-verified hairpins by transient transfection in 293T cells (if possible) or by transduction with lentiviruses as described in Janes_Virusprep.pdf. Coexpression of a V5-tagged expression construct from the ORFeome is highly recommended if there are concerns about antibody specificity.

shRNA cloning into pLKO.1 vectors

Janes Lab Protocols

Entered by Karin Holmberg, Kevin Janes

12/14/14

Buffer recipes

- **10x annealing buffer** Store at room temperature
 - 100 mM Tris-HCl (pH 7.5)
 - 1 M NaCl
 - 10 mM EDTA